

Millipore®

Preparation, Separation,
Filtration & Monitoring Products

Improved endotoxin removal using ecofriendly detergents

Intensified plasmid DNA capture

Juan Amor
DCVMN International 2024

MERCK

Agenda

Millipore®

Preparation, Separation,
Filtration & Monitoring Products



1- Plasmid Production overview

2- High productivity direct plasmid capture step

3- Endotoxin clearance strategy for plasmid capture

MERCK

Millipore®

Preparation, Separation,
Filtration & Monitoring Products

Plasmid Production Overview


01

MERCK

Background


Millipore®

What is plasmid and why do we care?

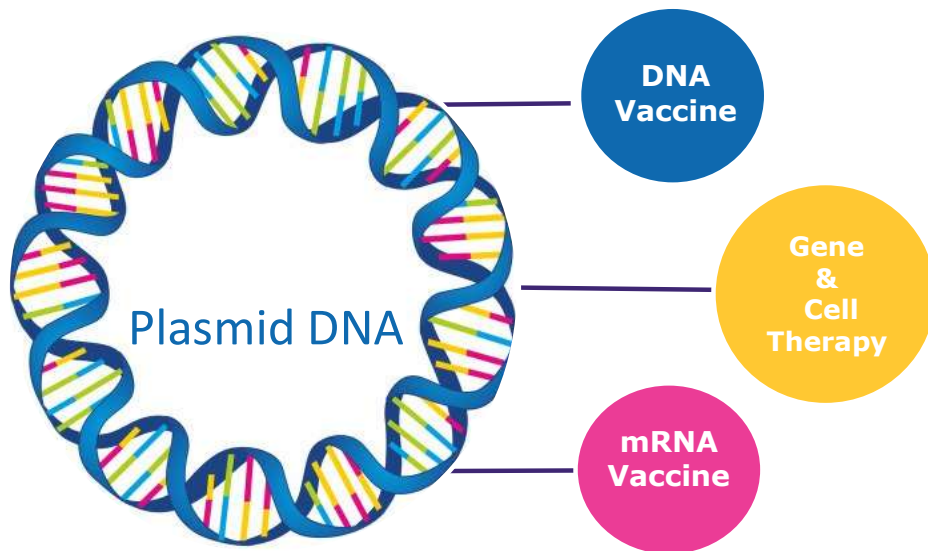

Characteristic

- **Extrachromosomal circular DNA**
- Commonly in **bacterial organisms**
- Sensitive to mechanical stress

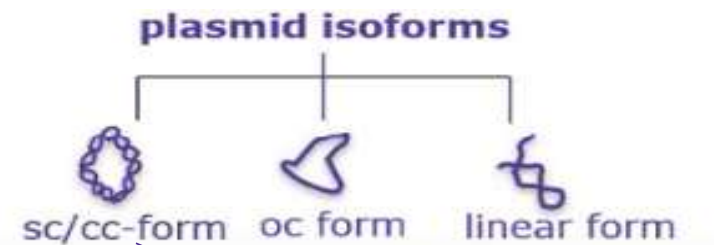

Therapeutically effective




Solutions to improve plasmid production are needed!

Process Overview


Plasmid Manufacturing

Millipore®

pDNA Manufacturing

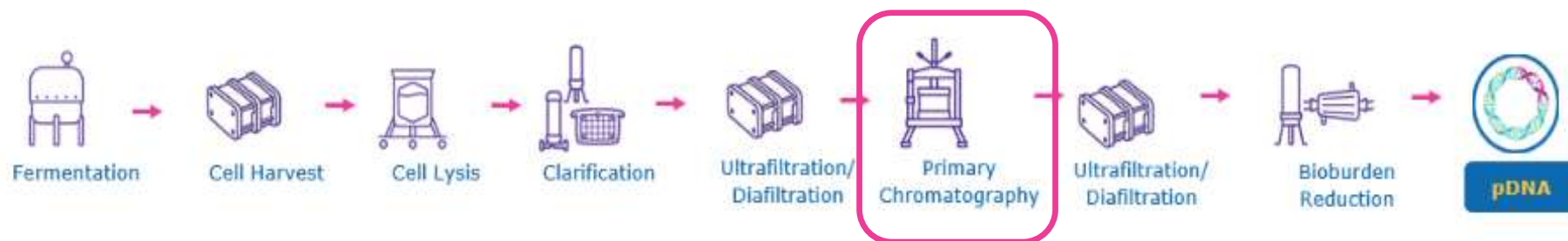

Chromatography is a current bottleneck

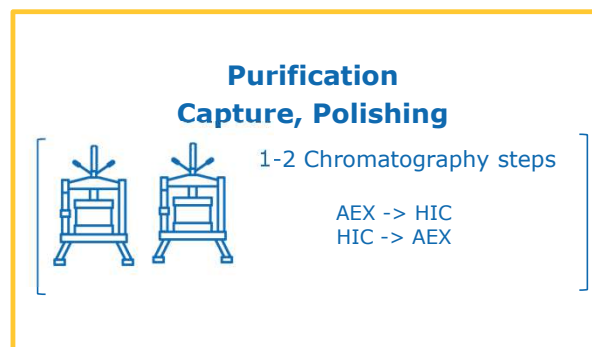

Millipore®

Preparation, Separation,
Filtration & Monitoring Products

High productivity direct plasmid capture step

02

MERCK



Natrix® Q Chromatography Membrane in Single-use Format

Improved performance enabling high-throughput capture

Millipore®

Material Properties

- Composite membrane, high mechanical strength
- Macroporous hydrogel, nominal pore size: 0.4 µm
- Large accessible surface for plasmids and other macromolecules

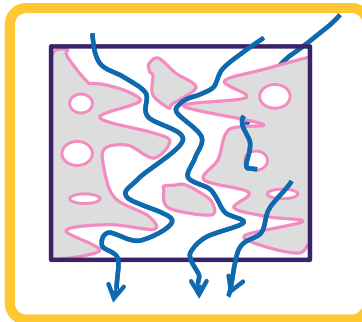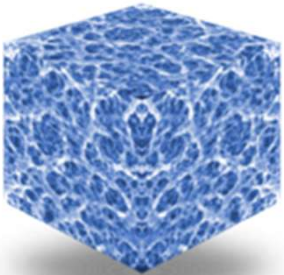

Plasmid DNA Case Study

Millipore®

AEX Capture Chromatography - Test Overview and Methods


Clarified lysate conditions:

- 6.5 kbp pDNA, 24 µg/mL titer. 1.5M K-acetate buffer, pH 5.3, 86.9 mS/cm
- Nucleic acid content: 3.8% pDNA, 96.2% RNA. Endotoxin content: 380,000 EU/mg pDNA



Step	Mobile Phase	Membrane Volumes	Flowrate
Equil	1M K-Acetate + 150 mM NaCl, pH 5.0 (75 mS/cm)	50 MV	10 MV/min
Load	Clarified, sterile filtered lysate pH 5.2	11 mg pDNA/mL membrane	10 MV/min
Wash	1M K-Acetate + 150 mM NaCl, pH 5.0 (75 mS/cm)	20 MV	10 MV/min
Elute	100 mM Tris, pH 9 + 1M NaCl	50 MV	5 MV/min
CIP	1M NaOH + 2 M NaCl	20 MV	10 MV/min

Impact of NaCl supplementation on RNA clearance

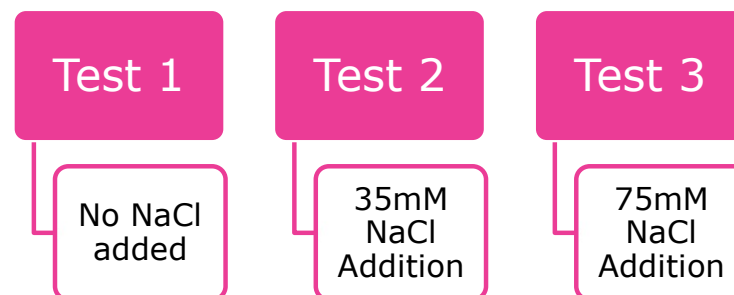

Capture pDNA while impurities (RNA) flowthrough

Analytics:

- DNA and RNA content assessed by HPLC (Tosoh DNA-NPR method)¹
- Endotoxin content assessed by Charles River Endosafe assay


¹ Urthaler 2005

Plasmid DNA Case Study

AEX Capture Chromatography - Results

Millipore®

Impact of Salt Supplementation on Capacity, Purity, Recovery

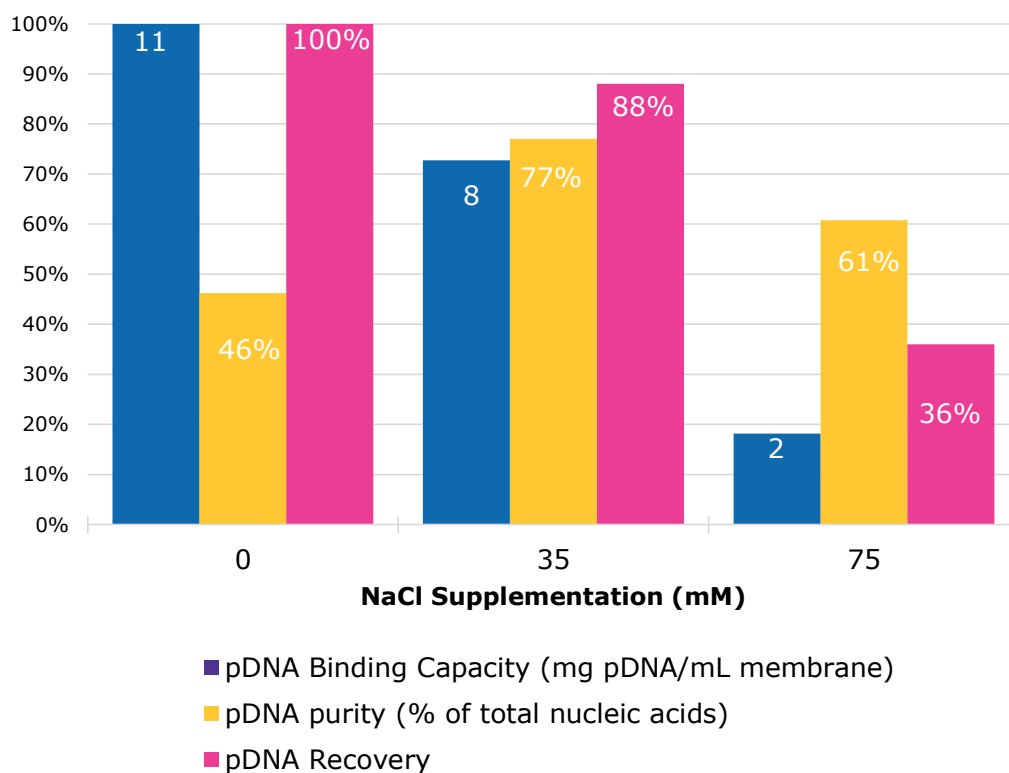

35 mM NaCl supplementation offers best balance of capacity, purity, recovery:

- Capacity = 8 mg pDNA/mL membrane
- Nucleic acid purity = 77% pDNA
- pDNA recovery = 88%
- Endotoxin content = 3,100 EU/mg pDNA



Can RNA and endotoxin purity be further improved with an alternative wash strategy?

Millipore®

Preparation, Separation,
Filtration & Monitoring Products

Endotoxin clearance strategy for plasmid capture

03

MERCK

Endotoxin

Molecule characteristics

Millipore®

- **Structure**

- Unique group of Lipopolysaccharides (LPS)
- Occur naturally in the cell wall of Gram-negative bacteria

- **Physico-Chemical Properties**

- Amphiphilic character: negatively charged core region <-> hydrophobic lipid A part
- Self-aggregating properties; formation of supramolecular structures (micells)

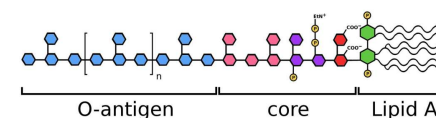


Fig 1. LPS molecular structure¹

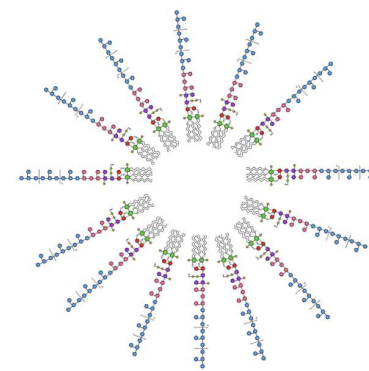


Fig 2. Micellar LPS structure

¹ Peri F, Pure Appl. Chem. 2012; 84(1): 97–106



Due to their pyrogenic effects, endotoxins are a critical impurity to remove from therapeutics.

Detergent Use in Bioprocessing

Solution to improve endotoxin clearance

- Some plasmid feeds may result in high endotoxin content in the AEX eluate
 - Dependent on cell culture characteristics, lysis conditions, AEX material, etc.
- **Triton™ X-100** has been used by the biopharmaceutical industry for effective endotoxin clearance
- **Triton™ X-100** has a broad range of uses:
 - Virus inactivation for blood-derived therapeutics, recombinant proteins
 - Cell lysis in viral vector production
 - Protein solubilization

However:

- As of January 2021, **TRITON™ X-100** has been prohibited in the EU unless authorization is granted
- This decision affects European manufacturers and downstream users



There is a need to identify eco-friendly TRITON™ X-100 alternatives

Detergents Use in Bioprocessing

Considerations when choosing a detergent

Millipore®

Process Requirements

- Effective endotoxin removal
- Does not adversely affect yield or quality of the product
- Low or no interference with product related analytical assays
- Does not interfere with subsequent purification
- Can be sufficiently removed from the product by the capture and subsequent DSP steps

Material Attributes

- Environmentally compatible
- High quality, purity
- Low toxicity potential

User Requirements

- Supply availability
- Suitability for different applications



Plasmid DNA Case Study



AEX Capture Chromatography- Wash Strategy

Control Wash

Step	Mobile Phase
Equilibration	1M K-Acetate + 150 mM NaCl, pH 5.0 (75 mS/cm)
Load	Clarified, sterile filtered lysate pH 5.2 + 35mM NaCl
Wash	1M K-Acetate + 150 mM NaCl, pH 5.0 (75 mS/cm)
Elute	100 mM Tris, pH 9 + 1M NaCl
CIP	1M NaOH + 2 M NaCl

Detergent Wash

Step	Mobile Phase
Equilibration	1M K-Acetate + 150 mM NaCl, pH 5.0 (75 mS/cm)
Load	Clarified, sterile filtered lysate pH 5.2 + 35mM NaCl
Wash	1M K-Acetate + 150 mM NaCl, pH 5.0 (75 mS/cm)
Detergent Wash	0.1M Tris, 10mM NaCl, + 0.5% detergent , pH 7.5
EDTA Wash	0.1M Tris, 10mM NaCl, + 2mM EDTA , pH 7.5
Elute w/ EDTA	100 mM Tris, 1M NaCl + 2mM EDTA , pH 9
CIP	1M NaOH + 2 M NaCl

Results

	Nucleic Acid Content	Endotoxin Content	Cycle Time
Feed Conditions	4% DNA, 96% RNA	380,000 EU/mg	N/A
Elution w/ Control wash (measured from eluate pool)	77% DNA, 23% RNA	3,100 EU/mg	55 min
Elution w/ Detergent wash (measured from eluate pool)	95% DNA, 5% RNA	500 EU/mg	65 min

Plasmid DNA Case Study

Millipore®

AEX Capture Chromatography- Membrane Vs Resin

100 L batch of clarified lysate, 3.6 g pDNA

Membrane Chromatography

Binding Capacity	8 g/L
Membrane Volume	0.46 L
Flow Rate	4.6 LPM
Step Time	1.1 hr
Cycles	1 cycle
Productivity	7.3 g pDNA/L/hr

Resin Chromatography

Binding Capacity	3 g/L
Resin Volume	1.18 L
Flow Rate	0.3 LPM
Step Time	9.9 hr
Cycles	1 cycle
Productivity	0.3 g pDNA/L/hr

Matrix® Q membrane is 24x more productive than resin



Conclusions

- 1 **Intensified plasmid capture solution: Membrane adsorber and optimal salt supplementation approach offers 3x plasmid binding capacity at 10x higher throughput.**
- 2 **Identified eco-friendly detergent for endotoxin removal that meet process, material, and user requirements**
- 3 **Identified eco-friendly detergent protocol for robust and high endotoxin clearance**
 - Integration into wash buffer preferred
 - Low residual detergent levels
 - No impact on plasmid binding capacity, yield, and purity
 - **Devirion™ C16** detergent: clearance >500x more than without detergent

Thank you

© 2024 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved. Merck, Millipore, SAFC, the vibrant M, Fractogel, Natrix, Tergitol, TWEEN, Triton and Deviron are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

